



Digital incubating/ refrigerating benchtop shakers.

Transparent lid offers great visibility and gas spring supports ensure hand-down handling of the samples by allowing immediate and easy access to your samples.



REMOTE CONTROL
with your MOBILE DEVICE.
see page 8-9



IST-3075 / IST-4075R

with optional dedicated platform and universal attachment

Optional accessories see page 95-97

- Universal platforms, Rubber mats
- Flask clamps, Funnel clamps, Test tube racks, Gateway

※ Above products design can be changed without prior notice.



Control display



Gas spring lid



Performance

- Temperature range:
 - Ambient +5°C to 80°C for IST-3075 / 4075.
 - Ambient -15°C (Min. 15) to 80°C for IST-3075R / 4075R.
- Competitive high shaking speed up to 500 rpm.
- Orbital shaking motion in 19mm diameter.
- Microprocessor PID control.
- Three point temperature calibration / Automatic tuning.
- Wide speed range even with heavy workload.
 - Equipped with a maintenance-free BLDC motor to generate smooth, quiet, uniform, and yet powerful shaking motion.
- Best effort run function intelligently manages its rpm to keep shaking even workload is out of its capacity.
 - Automatic shaking speed adjustment in case of excessive workload such as unbalanced load placement, unusual vibrations caused from unstable floor or external shock.
- Pleasant test environment by the smooth acceleration and deceleration control.
 - Smooth start and smooth stop function prevents chemical spills from flasks or test tubes.
- High-velocity fan ensures uniform temperature distribution and rapid thermal recovery after door open.
- Air-tight silicone door seal to provide excellent temperature uniformity.

Convenience

- Intuitive touch screen LCD controller with easy icons and logical menus.
- Easy-set digital timer for shaking operation.
(1 min. to 999 hr. 59 min.)
- Wide sample monitoring and convenient sample loading.
 - Convenient sample monitoring without affecting inside environment thanks to the three side transparent acrylic lid with bright LED lamps.
 - Gas spring lid supports providing smooth motion, soft opening and closing.
- Repetitive task can be easily performed as the platform stops where it starts.
 - Stopping the shaking platform always at the same position is highly beneficial for automated dosing or sampling process.
- Corrosion resistance stainless steel interior.
- Built-in RS-232 port and USB port for external control and data collection.

Safety

- Automatic run after power interruption.
- Temperature and shaking speed deviation alarm.
- Over-current protection, Stalled platform check.
- Mechanical over-temperature protection device.
- Environment-friendly CFC-free refrigerant. (for IST-3075R / 4075R)
- Audible and visible open door alarms.

Model		IST-3075		IST-3075R		IST-4075		IST-4075R	
Temperature	Range (°C / °F)	Amb. +5 to 80 / Amb. +9 to 176		Amb. -15 (Min. 15) to 80 / Amb. -27 (Min. 27) to 176		Amb. +5 to 80 / Amb. +9 to 176		Amb. -15 (Min. 15) to 80 / Amb. -27 (Min. 27) to 176	
	Fluctuation at 37°C in flask (±°C / °F)	0.1 / 0.18							
	Variation at 37°C in flask (±°C / °F)	0.5 / 0.90							
	Refrigerator (Hp)	-	1 / 6 HBP		-		1 / 6 HBP		
Shaking system	Motion type	Orbital							
	Amplitude size (mm / inch, dia.)	19.1 / 0.75							
	Speed range (RPM)	20 to 500							
	Accuracy	±1 % of set speed (>100rpm) / ±1 (≤100rpm)							
	Timer	1 min. to 999 hr 59 min.							
	Max. load (kg / lbs)	10 / 22.0 at 500 rpm 15 / 33.1 at 400 rpm				14 / 30.9 at 500 rpm 21 / 46.3 at 400 rpm			
Dimension	Volume (L / cu ft)	53 / 1.9				83 / 2.9			
	Platform (WxD) (mm / inch)	350×350 / 13.8×13.8				450×450 / 17.7×17.7			
	Interior (WxDxH) (mm / inch)	410×410×320 / 16.1×16.1×12.6				510×510×320 / 20.1×20.1×12.6			
	Exterior (WxDxH) (mm / inch)	440x785x510 / 17.3x30.9x20.1				540x 890 x510 / 21.3x35x20.1			
	Net weight (Kg / lbs)	65 / 143.3		73 / 160.9		81 / 178.6		90 / 198.4	
Electrical requirements (230V)		60Hz, 4A		50Hz, 4A		60Hz, 5.5A		50Hz, 5.5A	
Cat. No.		AAH23411K	AAH23422K	AAH23511K	AAH23522K	AAH23611K	AAH23622K	AAH23711K	AAH23722K
Electrical requirements (120V)		60Hz, 7.5A		60Hz, 10.1A		60Hz, 7.5A		60Hz, 10.1A	
Cat. No.		AAH23465U		AAH23565U		AAH23665U		AAH23765U	

※ FDA establishment registered company. FDA listed products.

Incubated Shakers (Floor Standing)

Large capacity up to five 6L flasks and heavy workload shaking with rapid heating and cooling performance.

Exclusive drive mechanism with ergonomic design delivers years of trouble-free service.



REMOTE CONTROL
with your **MOBILE DEVICE**
see page 8-9



ISF-7100R

Optional accessories see page 95-97

- Universal/Dedicated platforms
- Spring wire racks, Microplate holders, Gateway

※ Above products design can be changed without prior notice.



User-friendly design



Clear viewing window



Gas spring lid



Performance

- Temperature range:
 - Ambient +5°C to 80 °C for ISF-7100 / 7200
 - Ambient -20°C (Min. 4) to 80°C for ISF-7100R / 7200R
- Competitive high shaking speed up to 500 rpm.
 - from 30 to 500 rpm for ISF-7100(R)
 - from 30 to 300 rpm for ISF-7200(R)
- Orbital shaking motion in various diameter.
 - 25.4mm for ISF-7100(R)
 - 50.8mm for ISF-7200(R)
- Microprocessor PID control.
- Three point temperature calibration / Automatic tuning.
- Wide speed range even with heavy workload.
 - Ideal DD(Direct Drive) brushless motor and triple cam system support heavy load, quiet, maintenance free operation and deliver high torque at high speeds and smooth rotation.
- Best effort run function intelligently manages its rpm to keep shaking even workload is out of its capacity.
 - Automatic shaking speed adjustment in case of excessive workload such as unbalanced load placement, unusual vibrations caused from unstable floor or external shock.
- Pleasant test environment by the smooth acceleration and deceleration control.
 - Smooth start and smooth stop function prevents chemical spills from flasks or test tubes.
- High-velocity fan ensures uniform temperature distribution and rapid thermal recovery after door open.
- Air-tight silicone door seal with lockable door for safe storage of samples.

Convenience

- Intuitive touch screen LCD controller with easy icons and logical menus.

- Easy-set digital timer for shaking operation. (1 min. to 999 hr. 59 min.)
- Easy sample monitoring and convenient sample loading.
 - Clear observation of samples without affecting inner chamber's environment thanks to the transparent top viewing window with bright LED lamps.
 - The lid is inclined gently towards the front for easy observation of samples in the chamber.
 - Gas spring lid supports providing smooth motion, soft opening and closing.
- Repetitive task can be easily performed as the platform stops where it starts.
 - Stopping the shaking platform always at the same position is highly beneficial for automated dosing or sampling process.
- Corrosion resistance stainless steel interior.
- Easy-access drain system with a quick disconnect valve offers great convenience to clean up of spills.
- Built-in RS-232 port and USB port for external control and data collection.
- Sturdy foot caster, beneficial for easy mobility during installation or relocation.

Safety

- Automatic run after power interruption.
- Temperature and shaking speed deviation alarm.
- Over-current protection.
- Triple independent temperature monitor system.
 - Temperature deviation alarm (High/Low)
 - Electronic temperature limiter (High/Low)
 - Mechanical over-temperature limiter (High)
- Environment-friendly CFC-free refrigerant. (for ISF-7100R / 7200R)
- Audible and visible open door alarms.

Model		ISF-7100		ISF-7100R		ISF-7200		ISF-7200R					
Temperature	Range (°C / °F)	Amb. +5 to 80 / Amb. +9 to 176		Amb. -20 (Min. 4) to 80 / Amb. -36 (Min. 22) to 176		Amb. +5 to 80 / Amb. +9 to 176		Amb. -20 (Min. 4) to 80 / Amb. -36 (Min. 22) to 176					
	Fluctuation at 37°C in flask (±°C / °F)	0.1 / 0.18											
	Variation at 37°C in flask (±°C / °F)	0.5 / 0.90											
	Refrigerator (Hp)	-	1/6Hp		-		1/6Hp						
Shaking system	Motion type	Orbital											
	Amplitude size (mm / inch, dia.)	25.4 / 1				50.8 / 2							
	Speed range (rpm)	30 to 500				30 to 300							
	Accuracy	±1 % of set speed (>100rpm) / ±1 (≤100rpm)											
	Timer	1 min. to 999 hr 59 min.											
	Max. load (kg / lbs)	15 / 33.1 at 500 rpm 25 / 55.1 at 400 rpm 35 / 77.2 at 300 rpm				20 / 44.1 at 300 rpm 35 / 77.2 at 200 rpm							
Dimension	Volume (L / cu ft)	270 / 9.5											
	Platform (WxD) (mm / inch)	755×520 / 29.7×20.5											
	Interior (WxDxH) (mm / inch)	894×634×480 / 35.2×25×18.9											
	Exterior (WxDxH) (mm / inch)	1128×854×1035 / 44.4×33.6×40.8											
	Net weight (Kg / lbs)	277 / 610.7				287 / 632.7		277 / 610.7		287 / 632.7			
Electrical requirements (230V)		50 / 60Hz, 3.7A		60Hz, 5.6A		50Hz, 5.6A		50 / 60Hz, 3.7A		60Hz, 5.6A		50Hz, 5.6A	
Cat. No.		AAH23342K		AAH23542K		AAH23541K		AAH23442K		AAH23642K		AAH23641K	
Electrical requirements (120V)		60Hz, 7.3A		60Hz, 10A		60Hz, 7.3A		60Hz, 7.3A		60Hz, 10A		60Hz, 10A	
Cat. No.		AAH23375U		AAH23575U		AAH23475U		AAH23475U		AAH23675U		AAH23675U	

※ Temp. change may occur during automatic defrost. (for ISF-7100R/7200R)

※ FDA establishment registered company. FDA listed products.

Stackable up to three units for maximum utilization of lab spaces.

Multi-purpose incubated shakers feature a wide range of temperature control and large chamber volume.



REMOTE CONTROL
with your **MOBILE DEVICE**
see page 8-9



Advantages of Stackable Model

- Large shaking capacity chamber allows user to freely organize different size of samples.
- 198L chamber can accommodate six of 4L flasks and also thirty nine of 250ml flasks. You can make the capacity triple, if you stack three units.
- Maximize your space with our triple stackable incubated shaker.
- Highly recommended for users who want to maximize limited space. Have three units in your lab while taking up minimal space. In addition, high rpm is still guaranteed even when you stack up 3 units.
- Lift up opening door and slide out platform enables easy sample handling.
- Door is opened upward and stayed open for hands-free handling. Slide type universal platform let you access to all samples, even ones placed at the back.
- Optional stand is available and highly recommended for comfort use. Setting your incubated shaker at your height let you feel totally relaxed when you use it anytime. (optional)

ISS-7100R / ISS-7200

with optional accessories

Standard accessories

- Stacking kits, Retractable foot caster

Optional accessories

- Universal platforms, Flask clamps, Universal attachments
- Stands, Gateway

see page 95-97

※ Above products design can be changed without prior notice.

Control display



Lift up opening door





Performance

- Temperature range.
 - Ambient +10°C to 80°C for ISS-7100 / 7200.
 - Ambient -20°C (Min. 4) to 80°C for ISS-7100R / 7200R.
- Competitive high shaking speed up to 400 rpm. (Stackable up to 250rpm)
 - from 30 to 400 rpm (stackable : 30 to 250) for ISS-7100(R)
 - from 30 to 300 rpm (stackable : 30 to 200) for ISS-7200(R)
- Orbital shaking motion in various diameters.
 - 25.4mm for ISS-7100(R)
 - 50.8 mm for ISS-7200(R)
- Microprocessor PID control.
- Three point temperature calibration / Automatic tuning.
- Wide speed range even with heavy workload.
 - Ideal DD(Direct Drive) brushless motor and triple cam system support heavy load, quiet, maintenance free operation and deliver high torque at high speeds and smooth rotation.
- Best effort run function intelligently manages its rpm to keep shaking even workload is out of its capacity.
 - Automatic shaking speed adjustment in case of excessive workload such as unbalanced load placement, unusual vibrations caused from unstable floor or external shock.
- Pleasant test environment by the smooth acceleration and deceleration control.
 - Smooth start and smooth stop function prevents chemical spills from flasks or test tubes.
- High-velocity fan ensures uniform temperature distribution and rapid thermal recovery after door open.
- Air-tight silicone door seal to provide excellent temperature uniformity.

Convenience

- Intuitive touch screen LCD controller with easy icons and logical menus.
- Easy-set digital timer for shaking operation. (1 min. to 999 hr. 59 min.)
- Wide sample monitoring.
 - Clear observation of samples without affecting inner chamber's environment thanks to the transparent viewing window with bright LED lamps.
- Repetitive task can be easily performed as the platform stops where it starts.
 - Stopping the shaking platform always at the same position is highly beneficial for automated dosing or sampling process.
- Built-in electrical outlet with a safety cover inside the chamber.
- Corrosion resistance stainless steel interior.
- Easy-access drain system with a quick disconnect valve offers great convenience to clean up of spills.
- Built-in RS-232 port and USB port for external control and data collection.
- The unit comes with retractable foot caster, beneficial for easy mobility during installation or relocation.

Safety

- Automatic run after power interruption.
- Temperature and shaking speed deviation alarm.
- Over-current protection.
- Triple independent temperature monitor system.
 - Temperature deviation alarm (High/Low)
 - Electronic temperature limiter (High/Low)
 - Mechanical over-temperature limiter (High)
- Environment-friendly CFC-free refrigerant. (for ISS-7100R / 7200R)

Model		ISS-7100		ISS-7100R		ISS-7200		ISS-7200R	
Temperature	Range (°C / °F)	Amb. +10 to 80 / Amb. +18 to 176		Amb. -20 (Min. 4) to 80 / Amb. -36 (Min. 72) to 176		Amb. +10 to 80 / Amb. +18 to 176		Amb. -20 (Min. 4) to 80 / Amb. -36 (Min. 72) to 176	
	Fluctuation at 37°C in flask (±°C / °F)	0.1 / 0.18							
	Variation at 37°C in flask (±°C / °F)	0.5 / 0.90							
	Refrigerator (Hp)	-	1/6Hp		-		1/6Hp		
Shaking system	Motion type	Orbital							
	Amplitude size (mm / inch, dia.)	25.4 / 1				50.8 / 2			
	Speed range (rpm)	30 to 400 (stackable ² : 30 to 250)				30 to 300 (stackable ² : 30 to 200)			
	Accuracy	±1% of set speed (>100rpm) / ±1 (≤100rpm)							
	Timer	1 min. to 999 hr 59 min.							
	Max. Load (kg / lbs)	20 / 44.1 at 400 rpm 15 / 33.1 at 250 rpm (stacked) 20 / 33.1 at 200 rpm (stacked)				20 / 44.1 at 300 rpm 15 / 33.1 at 200 rpm (stacked) 20 / 44.1 at 150 rpm (stacked)			
Dimension	Volume (L / cu ft)	198 / 7.0							
	Platform (WxD, mm / inch)	755x520 / 29.7x20.5							
	Interior (WxDxH, mm / inch)	809x624x393 / 31.9x24.6x15.5							
	Exterior ¹⁾ (WxDxH, mm / inch)	1310x996x755 / 51.6x39.2x29.7							
	Net weight (kg / lbs)	365 / 804.7		375 / 826.7		365 / 804.7		375 / 826.7	
Electrical requirements (230V)		60Hz, 3.7A 50Hz, 3.7A		60Hz, 5.6A 50Hz, 5.6A		60Hz, 3.7A 50Hz, 3.7A		60Hz, 5.6A 50Hz, 5.6A	
Cat. No.		AAH238131K AAH238132K		AAH238231K AAH238232K		AAH239131K AAH239132K		AAH239231K AAH239232K	
Electrical requirements (120V)		60Hz, 7.1A		60Hz, 10A		60Hz, 7.1A		60Hz, 10A	
Cat. No.		AAH238115U		AAH238215U		AAH239115U		AAH239215U	

※ Temp. change may occur during automatic defrost. (for ISS-7100R/7200R) ※ FDA establishment registered company. FDA listed products.

1) Main body dimension with retractable foot caster.

2) Stacked with stand or products.

Specially designed to be double-stacked on the floor or located on the lab benches for space saving.

Multi-purpose incubated shakers feature a wide range of temperature control and large chamber volume.



REMOTE CONTROL
with your **MOBILE DEVICE**
see page 8-9



ISS-4075R / ISS-3075

with optional accessories

Standard accessories

- Wire shelves (2ea)

Optional accessories

see page 95-97

- Universal platforms, Universal attachments, Flask clamps
- Retractable foot caster, Stacking kits for stackable chambers, Gateway

※ Above products design can be changed without prior notice.

Control display





Performance

- Temperature range:
 - Ambient +5°C to 80°C for ISS-3075 / 4075.
 - Ambient -20°C (Min. 4) to 80°C for ISS-3075R / 4075R.
- Competitive high shaking speed up to 500 rpm.
(Stackable up to 250rpm)
- Orbital shaking motion in 19.1mm diameter.
- Microprocessor PID control.
- Three point temperature calibration / Automatic tuning.
- Wide speed range even with heavy workload.
 - Equipped with a maintenance-free BLDC motor to generate smooth, quiet, uniform, and yet powerful shaking motion.
- Best effort run function intelligently manages its rpm to keep shaking even workload is out of its capacity.
 - Automatic shaking speed adjustment in case of excessive workload such as unbalanced load placement, unusual vibrations caused from unstable floor or external shock.
- Pleasant test environment by the smooth acceleration and deceleration control.
 - Smooth start and smooth stop function prevents chemical spills from flasks or test tubes.
- High-velocity fan ensures uniform temperature distribution and rapid thermal recovery after door open.
- Air-tight silicone door seal to provide excellent temperature uniformity.

Convenience

- Intuitive touch screen LCD controller with easy icons and logical menus.
- Easy-set digital timer for shaking operation. (1 min. to 999 hr. 59 min.)

- Wide sample monitoring.
 - Clear observation of samples without affecting inner chamber's environment thanks to the transparent viewing window with bright LED lamps.
- Repetitive task can be easily performed as the platform stops where it starts.
 - Stopping the shaking platform always at the same position is highly beneficial for automated dosing or sampling process.
- Two adjustable-height shelves provided allow static incubation or refrigeration, which increase storage capacity.
- Built-in electrical outlet with a safety cover inside the chamber.
- Corrosion resistance stainless steel interior.
- Easy-access drain system with a quick disconnect valve offers great convenience to clean up of spills.
- Built-in RS-232 port and USB port for external control and data collection.
- Retractable foot caster, beneficial for easy mobility during installation or relocation. (optional)
- Chambers stackable, up to two levels.

Safety


- Automatic run after power interruption.
- Temperature and shaking speed deviation alarm.
- Over-current protection, Stalled platform check.
- Triple independent temperature monitor system.
 - Temperature deviation alarm (High/Low)
 - Electronic temperature limiter (High/Low)
 - Mechanical over-temperature limiter (High)
- Environment-friendly CFC-free refrigerant. (for ISS-3075R / 4075R)
- Audible and visible open door alarms.

Model		ISS-3075	ISS-3075R	ISS-4075	ISS-4075R
Temperature	Range (°C / °F)	Amb. +5 to 80 / Amb. +9 to 176	Amb. -20 (Min. 4) to 80 / Amb. -36 (Min. 7.2) to 176	Amb. +5 to 80 / Amb. +9 to 176	Amb. -20 (Min. 4) to 80 / Amb. -36 (Min. 7.2) to 176
	Fluctuation at 37°C in flask (±°C / °F)	0.1 / 0.18			
	Variation at 37°C in flask (±°C / °F)	0.5 / 0.90			
	Refrigerator (Hp)	-	1/6Hp	-	1/6Hp
Shaking system	Motion type	Orbital			
	Amplitude size (mm / inch, dia.)	19.1 / 0.75			
	Speed range (rpm)	20 to 500 (stackable : 20 to 250)			
	Accuracy	±1% of set speed (>100rpm) / ±1 (≤100rpm)			
	Timer	1 min. to 999 hr 59 min.			
	Max. Load (kg / lbs)	10 / 22.0 at 500 rpm 15 / 33.1 at 400 rpm 15 / 33.1 at 250 rpm (stacked)		10 / 22.0 at 500 rpm 21 / 46.3 at 400 rpm 21 / 46.3 at 250 rpm (stacked)	
Dimension	Volume (L / cu ft)	80 / 2.8		150 / 5.3	
	Platform (WxD, mm / inch)	350×350 / 13.8×13.8		450×450 / 17.7×17.7	
	Interior (WxDxH, mm / inch)	440×440×418 / 17.3×17.3×16.5		540×540×518 / 21.3×21.3×20.4	
	Exterior (WxDxH, mm / inch)	578×815×854 / 22.8×32.1×33.6		678×895×934 / 26.7×35.2×36.8	
	No. of wire shelf (standard / max.)	2 / 8		2 / 11	
Electrical requirements (230V)		60Hz, 6.2A 50Hz, 6.2A		60Hz, 6.2A 50Hz, 6.2A	
Cat. No.		AAH23431K AAH23432K		AAH23631K AAH23632K	
Electrical requirements (120V)		60Hz, 9.5A		60Hz, 9.5A	
Cat. No.		AAH23415U		AAH23615U	
		AAH23515U		AAH23715U	

※ Temp. change may occur during automatic defrost. (for ISS-3075R/4075R)

※ FDA establishment registered company. FDA listed products.

for Incubated Shakers

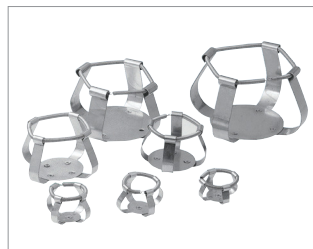

Flask clamps



Plastic clamps (plastic)



Funnel clamps



Dedicated platforms

Universal platforms / Flask clamps / Funnel clamps

Universal platforms provide flexibility for mixing assortments of test tubes and glassware of varying sizes on a single platform. Various clamps for Erlenmeyer flasks and separatory funnels are available.

Description		Model	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
		Platform size	350x350mm / 13.8x13.8inch	450x450mm / 17.7x17.7inch	755x520mm / 29.7x20.5inch	755x520mm / 29.7x20.5inch	350x350mm / 13.8x13.8inch	450x450mm / 17.7x17.7inch
Universal platform			AAA23501-V3	AAA23502-V3	AAA23503-V3	AAA23503-V4	AAA23501-V3	AAA23502-V3
Flask clamps (Stainless steel)	Erlenmeyer flask size	Max. mountable flask clamps						
	for 50 ml	AAA23550	36	49	88	88	36	49
	for 100 ml	AAA23551	24	39	68	68	24	39
	for 250 ml	AAA23552	14	20	39	39	14	24
	for 300 ml	AAA23556	14	20	39	39	14	24
	for 500 ml	AAA23553	9	16	28	28	9	16
	for 1L	AAA23554	6	9	18	18	6	9
	for 2L	AAA23555	4	5	11	11	4	6
	for 2.8L	AAA23557	2	4	8	8	2	5
	for 4L	AAA23558	-	-	6	6	-	4
	for 6L	AAA23559	-	-	5	-	-	2
Flask clamps (plastic)	Erlenmeyer flask size	Max. mountable flask clamps						
	for 50 ml	AAA30570	36	49	88	88	36	49
	for 100 ~ 125 ml	AAA30571	19	29	48	48	19	29
	for 200 ml	AAA30572	13	21	35	33	13	21
	for 250 ml	AAA30573	13	19	35	33	13	19
	for 300 ml	AAA30574	12	19	35	33	12	19
	for 500 ml	AAA30575	9	12	24	24	9	12
	for 1L	AAA30576	4	9	12	12	5	9
	for 2L	AAA30577	2	4	8	8	4	5
Funnel clamps	separatory funnel size	Max. mountable funnel clamps						
	for 250 ml	AAA23562	4	6	15	15	4	6
	for 500 ml	AAA23563	2	4	10	10	3	4
	for 1L	AAA23564	-	3	7	7	-	3
	for 2L	AAA23565	-	-	5	5	-	-

※ Above accessories need to be mounted on a universal platform which is sold separately.

Dedicated platforms

When using a single flask size only, the dedicated platforms with pre-installed clamps provide maximum capacity. 250mL and 500mL clamps are standard but other sizes can be available.

Description	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
Universal platform & 250mL flask clamps	AAA23611-V1	AAA23623-V1	AAA23634	AAA23637	AAA23611-V1	AAA23621-V1
No. of flask clamps	14	20	39	39	14	20
Universal platform & 500mL flask clamps	AAA23612-V1	AAA23622-V1	AAA23635	AAA23636	AAA23612-V1	AAA23622-V1
No. of flask clamps	9	16	28	28	9	16


* A dedicated platform is sold with setting each max. mountable amount of flask clamps on.

Accessories & Options


for Incubated Shakers



Test tube racks

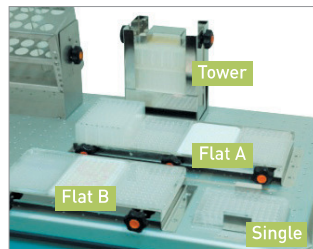

Microplate holders
(universal platform is optional)

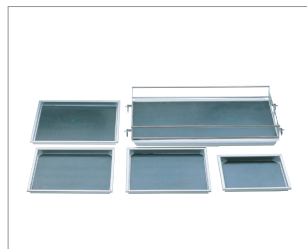

Rubber mats



Universal attachments

Test tube racks

Easy to release and mount racks made of durable stainless steel with adjustable angle and tube height support. Test tube racks are used with an optional universal platform.

Description	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
Universal platform	AAA23501-V3	AAA23502-V3	AAA23053-V3	AAA23503-V4	AAA31501-V3	AAA23502-V3
Test tube size	Mountable capacity of rack*					
for 86 test tubes, ø8mm / ø0.3"	2	4	7	7	2	4
for 86 test tubes, ø10mm / ø0.4"						
for 58 test tubes, ø12mm / ø0.5"						
for 58 test tubes, ø14mm / ø0.6"						
for 32 test tubes, ø17mm / ø0.7"						
for 19 test tubes, ø25mm / ø1.0"						
for 10 test tubes, ø30mm / ø1.2"						
for 10 test tubes, ø35mm / ø1.4"						
for 4 test tubes, ø50mm / ø1.9"						

* No declination

※ We can manufacture other size of test tube racks to meet your specific requirements. Contact us for details.

Microplate holders

Available in three styles: tower type, flat type (small and large), and single type. Tower type holder is for 10 standard or deep well plates. Large flat holder is for 4 standard or deep well plates. Small flat holder is for 3 standard or deep well plates. Single type holder is for 1 standard or deep well plate. (standard well plates' height: 14.7mm)

Description	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
Universal platform	AAA23501-V3	AAA23502-V3	AAA23053-V3	AAA23503-V4	AAA31501-V3	AAA23502-V3
Microplate holder type	Max. mountable microplate holders					
Single	AAA23654	8	15	25	8	15
Tower	AAA23651	5	8	21	6	10
Flat A(large)	AAA23652	-	2	4	-	2
Flat B(small)	AAA23653	2	2	6	2	3

* Microplate holders need to be mounted on an universal platform which is sold separately.

Rubber mats

Designed for low-speed applications, non-skid rubber mats allow quick addition or removal of flasks, plates or tubes.

Description	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
Cat. No.	AAA23531	AAA23532	AAA23534	AAA23535	AAA31531	AAA31532


Universal attachments (utility carriers)

Several adjustable clamping bars for quickly securing a variety of glassware or tubes on non-skid rubber mats. Suitable for flat-bottomed flasks of various sizes.

Description	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
Cat. No.	AAA23511-V1	AAA23512-V1	AAA23505	AAA23513	AAA23511-V1	AAA23512-V1



Spring wire racks
(two tier rack with flexible crossed spring)

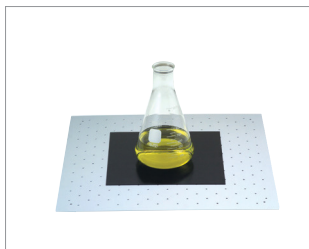

Lab sticker
(universal platform is optional)



Stand (stand 1 / stand 2)



Retractable foot caster

Spring wire racks

Ideal for holding a wide variety of tubes, flasks, and other glassware. All components are made of stainless steel.

Description	IST-3075(R)	IST-4075(R)	ISF-7100(R) ISF-7200(R)	ISS-7100(R) ISS-7200(R)	ISS-3075(R)	ISS-4075(R)
Cat. No.	AAA23521-V1	AAA23522-V2	AAA23514	AAA23526	AAA31521-V1	AAA31522-V1
Erlenmeyer flask size	Max. inserting flasks					
for 50 mL	16	25	45	45	16	25
for 100 mL	9	16	28	28	9	16
for 250 mL	4	9	18	18	4	9
for 300 mL	4	9	18	18	4	9
for 500 mL	4	5	15	15	4	5
for 1L	2	4	10	10	2	4
for 2L	1	2	6	6	1	2
for 2.8L	1	2	4	4	1	2

Lab sticker

An alternative method for securing vessels to a shaker platform. Ideal for containers such as bottles and dishes that cannot be held on the universal platform with conventional flask clamps. For safe operation, gradual RPM increase is recommended.

Cat. No.	Description (WxDxH) (mm / inch)
AAA30551	200×200×5 / 7.9×7.9×0.2

Stacking kits

Units are stackable using exclusive stacking kits for space saving.

Cat. No.	Description	Suitable for
AAA23573	Stacking kits for stackable chambers	ISS-7100(R), ISS-7200(R)
AAA23811	Stacking kits for stackable chambers	ISS-3075(R)
AAA23812	Stacking kits for stackable chambers	ISS-4075(R)

Stands

Perfectly designed to let users feel totally relaxed by leveling the height.

Cat. No.	Description	Suitable for
AAA23571	Stand 1 with Stacking kits (600mm / 23.6" height for stacking 1 unit)	ISS-7100(R), ISS-7200(R)
AAA23572	Stand 2 with Stacking kits (300mm / 11.8" height for stacking 2 units)	

Retractable foot caster

Highly beneficial for easy mobility during installation or relocation.

Cat. No.	Description	Suitable for
AAA23801	Retractable foot caster (700×733×145mm / 27.6×28.9×5.7")	ISS-3075(R)
AAA23802	Retractable foot caster (800×813×225mm / 31.5×32×8.9")	ISS-4075(R)

Gateway

Establish Jeio Tech Network by Gateway. No computer use is required at all. Remote monitoring of operating status through a mobile device (Android only) is simply available when you place a Gateway in your laboratory. Maximum 4 units can be connected to Gateway by RS-232 cable.

Cat. No.	Description	Suitable for
AAAQ1011	Gateway (2.4GHz, 100~240V, 50/60Hz)	IL3 models IST, ISF, ISS models



Gateway



Remote control through a mobile device
(Android only)